Installation

S19-949

**On-Site® Gravity-Fed Eye
Wash Refill Kit**

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Installation

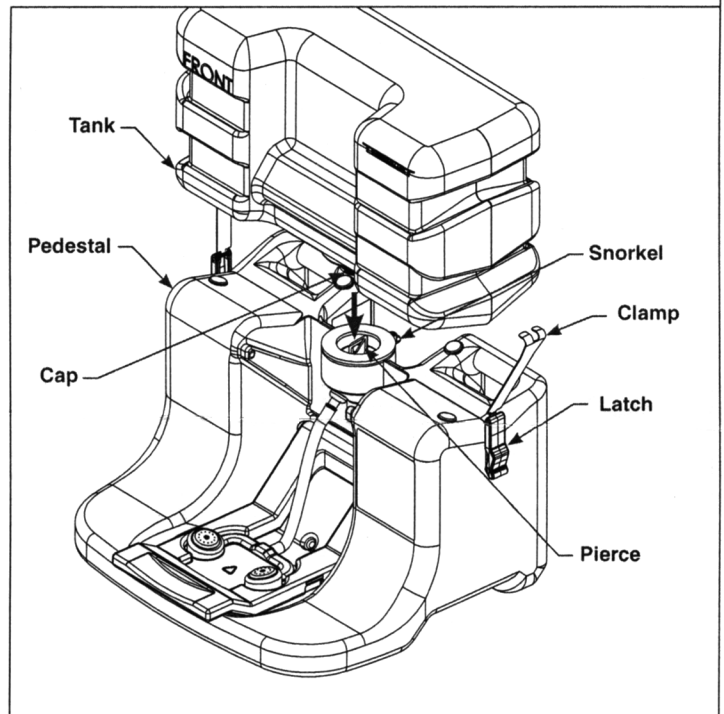
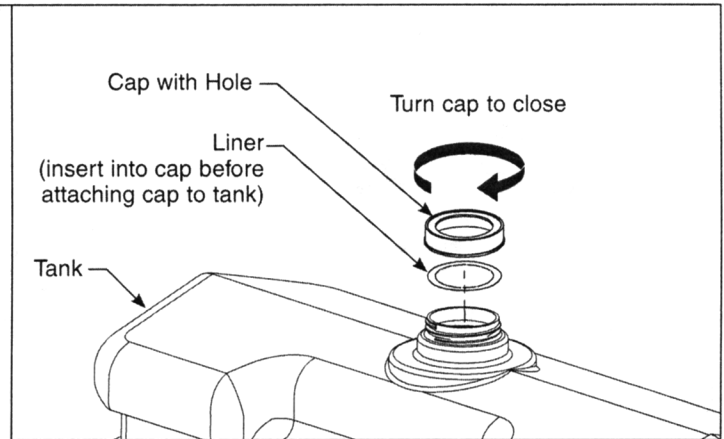
Fill and mount tank

1. Pour clean, potable water into the tank; continue pouring until the tank is $\frac{1}{3}$ to $\frac{1}{2}$ full.
2. Add water preservative (not included) to the eyewash tank.
3. Add clean, potable water to the tank until the water level reaches the top of the tank.
4. Insert one liner into the cap with hole and attach the cap to the top of the eye wash tank (make sure the cap is tightly closed).

NOTE FOR TANK INSTALLATION: To ensure proper installation, the tank should be placed on the pedestal from the front of the unit.

5. Invert the tank and place it on top of the pedestal. The pierce on the pedestal will puncture a hole in the tank cap liner and allow the solution to flow.
6. Raise the arm to snap fit in an upright position, then secure the tank to the pedestal using the side tank clamps. Make sure the latches are pulled down tight to the pedestal.

NOTE: After the tank is mounted to the pedestal, minor spillage could occur. Allow the fluid to stabilize completely before inspecting the unit for leaks.

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